



Software Update

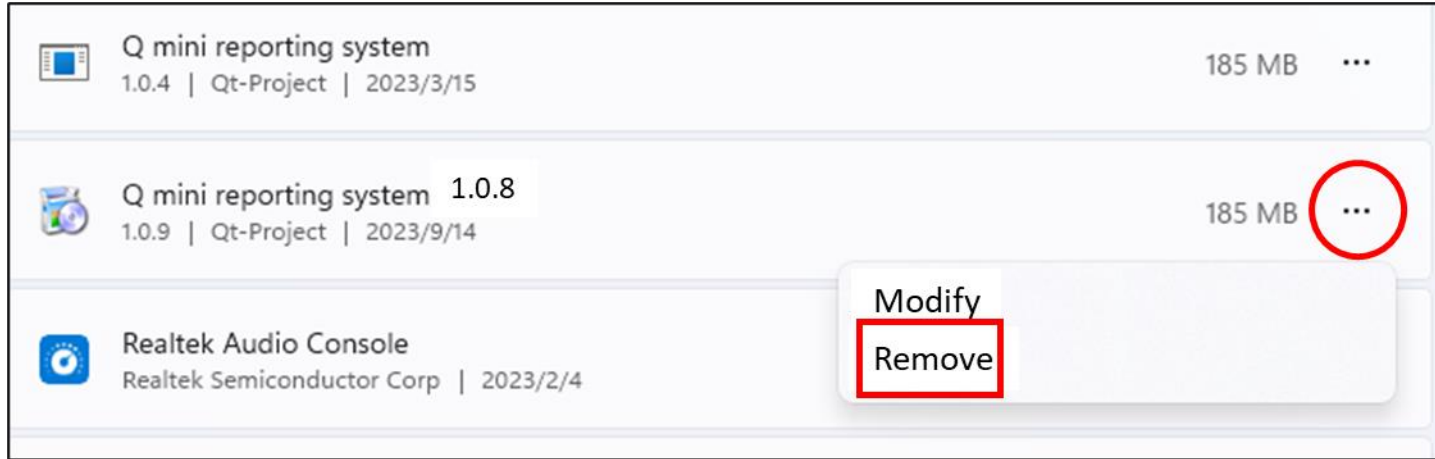

Step 1 :


Remove the old version software from the computer.

(If the software is not yet installed, simply download the new version file for installation.)

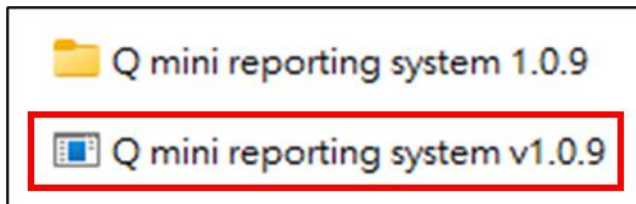


Step 2 :

Select the new software version for installation.

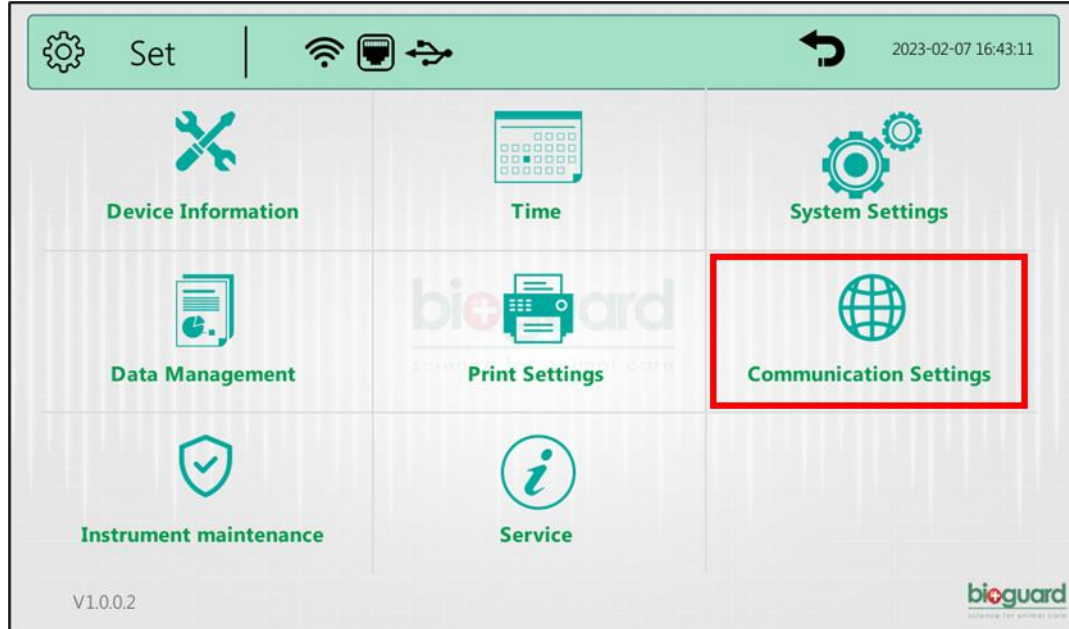

Step 3 :

Choose "Communication Settings"

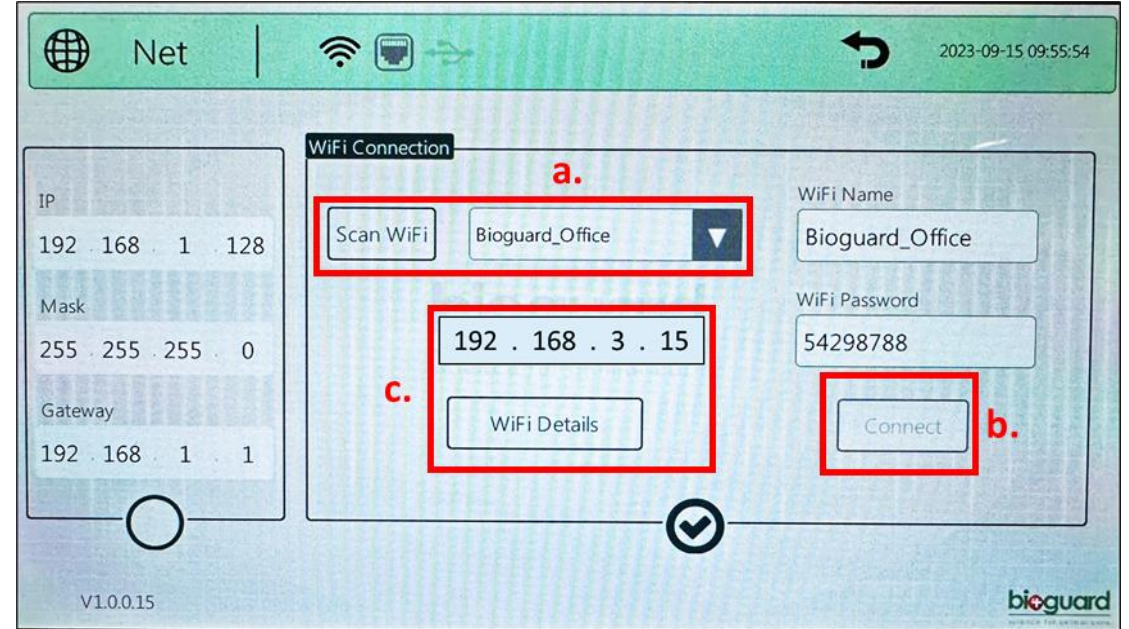

Step 4 :

a. Scan WiFi (Must be connected to the same network as the computer)


b. Click "Connect"

c. Click "WiFi Details" then show WiFi IP




Step 5 :

Start the newly installed software on the desktop.

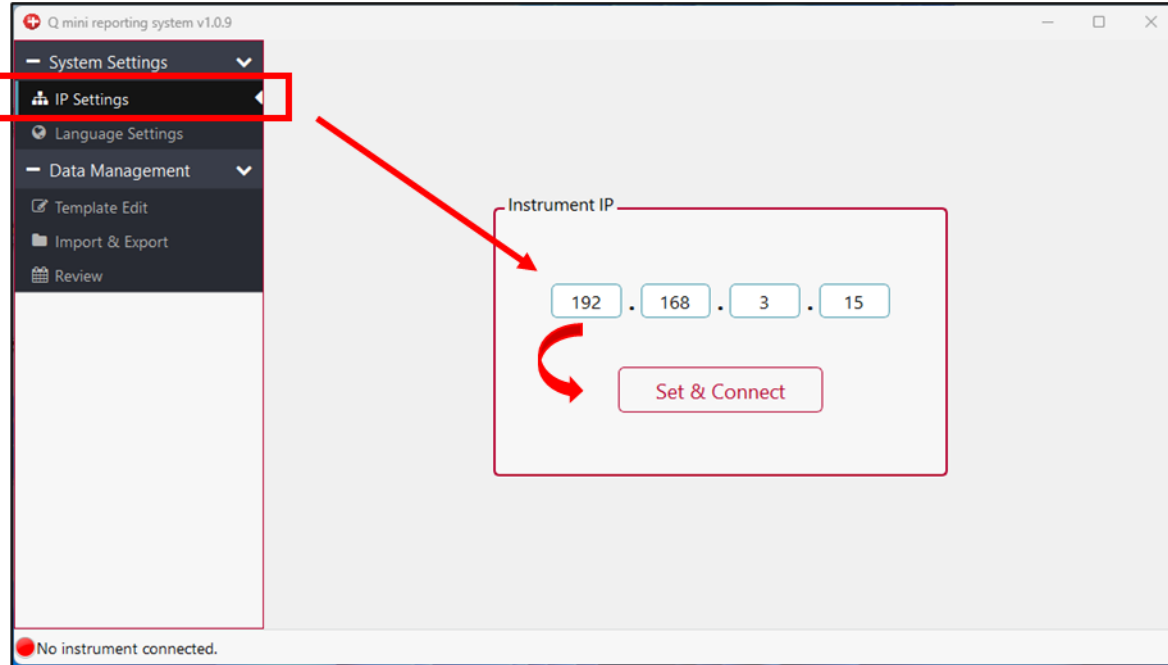

Step 6 :

Enter the IP information of the instrument.

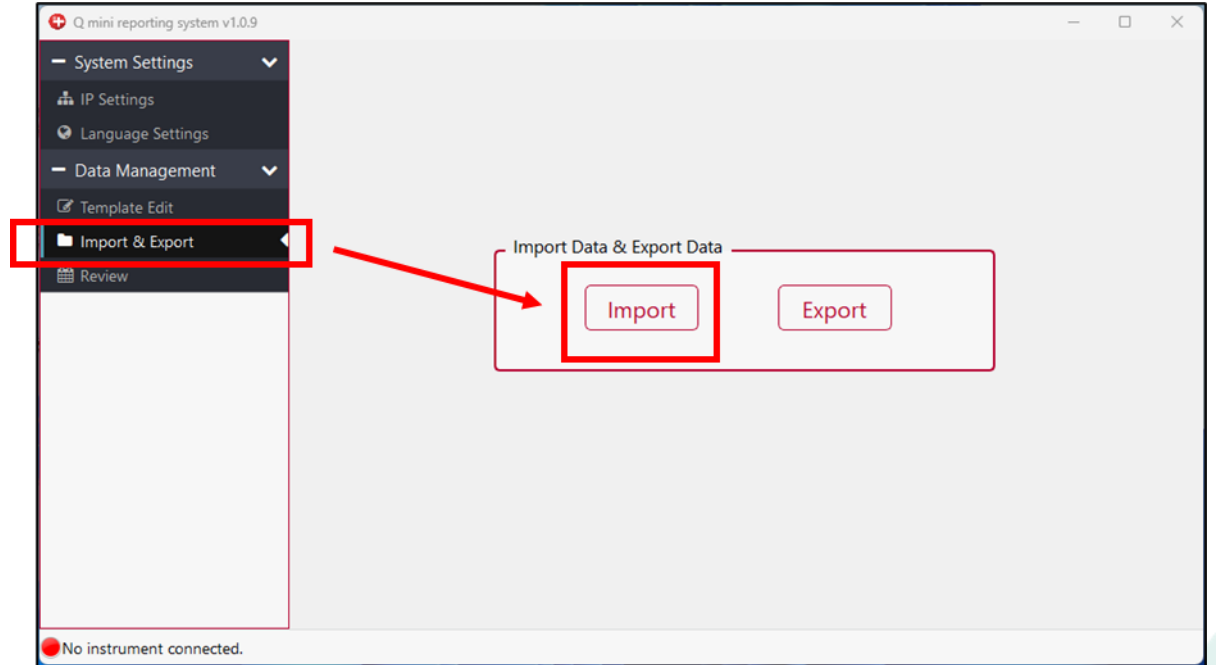

Step 7 :

Import old data from the device.